Product Use Limitation & Warranty

This product is intended to be used for life science research only. It has not been approved for drug or diagnostic purpose. YEASTERN's products should not be resold, modified for resale, or used to manufacture commercial products without written approval by YEASTERN. YEASTERN guarantees the performance of all products in the manner described in our protocol. The purchaser must determine the suitability of the product for its particular use. Should any product fail to perform satisfactorily due to any reason other than misuse, YEASTERN will replace it free of charge.

Ver. L0307

No part of these protocols may be reproduced in any form or by any mean, transmitted, or translated into a machine language without the permission of YEASTERN BIOTECH CO., LTD.

> Address: 6F-3, 23 Lane 169, Kang Ning St. Shijr, Taipei, 22180 Taiwan.

Email: veastern@veastern.com.tw





Copyright® 2011 All rights reserved. Yeastern Biotech Co., Ltd

Cat. No. FYT501-50R FYT501-100R

Concentration: 200 U/ul Storage: - 20 °C

Description

Deoxy* HiSpec Reverse Transcriptase (RT) is genetically engineered by introducing point mutations to MMLV RT that increase half-life, reduce RNase activity and increase thermal stability. Those designed mutations lead to increased specificity of Deoxy* HiSpec RT and the highest cDNA yield of all RTs. It is ideal for RT-PCR of a specific gene or generating cDNA from total or poly (A)* RNA samples. It synthesizes a complementary DNA strand from total RNA, mRNA, or an RNA:DNA hybrid.

Content

- Deoxy⁺ HiSpec Reverse Transcriptase
- 2X Deoxy⁺ RT premix:
 100 mM Tris HCl pH 8 3

100 mM Tris-HCl pH 8.3 \cdot 150 mM KCl \cdot 6 mM MgCl $_{\! 2}$ \cdot 20 mM DTT \cdot 1 mM dNTPs

Unit Definition

One unit incorporates 1 nmole of dTTP into acid precipitable material in 10 min at 37° C using poly(A)-oligo(dT) as template primer.


Standard Protocol for First-Strand cDNA Synthesis

Add the following components to the microtubes on ice

Oligo (dT) primer		50 pmole	
or Random primer		50 pmole	=
or Gene specific p	rimer	2 pmole	
2× Deoxy+ RT pr	emix	10 µl	
Template RNA	(total RNA ≤ 5µ	ug or mRNA ≤ 1μg)	
ddH₂O		variable	
	Oligo (dT) primer or Random primer or Gene specific p 2× Deoxy ⁺ RT pr Template RNA	Oligo (dT) primer or Random primer or Gene specific primer ≥ X Deoxy* RT premix Template RNA (total RNA ≤ 5)	or Random primer 50 pmole or Gene specific primer 2 pmole 2 × Deoxy+ RT premix 10 μ Template RNA (total RNA ≤ 5μg or mRNA ≤ 1μg)

Total volume 18 µl



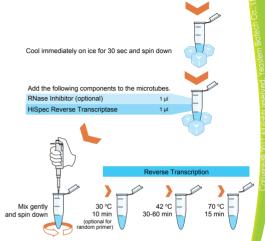

PCR (Recommended)

Use only 2 µl of the first-strand reaction for PCR.

. Add the following components to a PCR tu	be.
10× PCR Buffer 5	μl
10 mM dNTPs Mixture 1	μl
10 µM Forward primer 1	μl
10 µM Reverse primer 1	μl
5 U/µl Taq DNA polymerase 1	μl
The first-strand reactant 2	μΙ
ddH ₂ O to 50	μl

- 2. Mix gently and spin down.
- 3. Perform 20 to 40 cycles of PCR.