



Genomic DNA Extraction Mini Kit (Blood and Urine)

Product Use Limitation & Warranty

This product is intended to be used for life science research only. It has not been approved for drug or diagnostic purpose. YEASTERN's products should not be resold, modified for resale, or used to manufacture commercial products without written approval by YEASTERN. YEASTERN guarantees the performance of all products in the manner described in our protocol. The purchaser must determine the suitability of the product for its particular use. Should any product fail to perform satisfactorily due to any reason other than misuse, YEASTERN will replace it free of charge.

Ver. Q0306

No part of these protocols may be reproduced in any form or by any mean, transmitted, or translated into a machine language without the permission of YEASTERN BIOTECH CO., LTD.

Address: 6F-3, 23 Lane 169, Kang Ning St., Shijr, Taipei, 22180 Taiwan.
Tel: +886-2-2695-3922 **Fax:** +886-2-2695-3979
Email: yeastern@yeastern.com **Website:** www.yeastern.com

Copyright© 2017 All rights reserved. Yeastern Biotech Co., Ltd.
Copyright© 2017 All rights reserved. Yeastern Biotech Co., Ltd.

Cat. No.
SYG109-002
FYG109-100

Genomic DNA Extraction Mini Kit (Blood and Urine)

(SYG109-002/FYG109-100) Store at RT/4 °C Ver. Q0306

Sample : 300 µl (Whole Blood) /200 µl Buffy Coat /Urine Yield : Up to 50 µg

Contents

	SYG109-002	FYG109-100
B1 Buffer	4 ml	100 ml
B2 Buffer	1.5 ml	35 ml
B3 Buffer	0.5 ml	12 ml
BC Buffer	2 ml	45 ml
W1 Buffer	2 ml	45 ml
W2 Buffer*	0.3 ml x2	15 ml
Elution Buffer	1 ml	10 ml
BZ Column	4 pcs	100 pcs
Collection Tube	4 pcs	100 pcs
User Manual	1	1

Buffer Preparation

- Add ethanol (96-100%) to the W2 Buffer prior to first use

	SYG109-002	FYG109-100
W2 Buffer	300 ul x2	15 ml
ethanol (96 ~ 100%)	1.2 ml x2	60 ml

Important Notes

1. Buffers provided in this system contain irritants. Wear gloves and lab coat when handling these buffers.
2. Add ethanol (96- 100 %) to W2 Buffer when first open.
3. Prepare dry bath or water bath before the operation.
4. Resolve any precipitate by warming at 37 °C.

Purification Protocols

I. For Blood :

Step 1. Sample preparation

- a. Transfer up to 300 μ l of whole Blood or 200 μ l Buffy Coat to a microcentrifuge tube (not provided). (If sample volume is less than 200 μ l, adjust the volume to 200 μ l with PBS).

Step 2. RBC Lysis

- a. Add 900 μ l B1 Buffer to the tube and mix by inverting the tube. Incubate the mixture at room temperature for 10 minutes. During incubation, invert the tube every 5 minutes.

Step 3. Cell Lysis

- a. Centrifuge 5 mins at 4000 x g. Remove the supernatant completely.
- b. Resuspend the cell with 50 μ l B1 Buffer by pipetting the pellet.
- c. Add 300 μ l B2 Buffer to the sample and mix thoroughly by vortexing.
- d. Incubate at 60 °C for 10 mins until the sample lysate is clear. Invert the tube every 3 minutes during incubation.
- e. Preheat required Elution Buffer (50-90 μ l per sample) in a 60 °C water bath.
- f. Centrifuge 5 mins at 3000 x g. Remove the supernatant completely.

Optional Step

If RNA-free genomic DNA is required, perform this optional step.

- a. Add 5 μ l of RNase A (10 mg/ ml, not provided) to sample lysate and mix by vortexing.
- b. Incubate at room temperature for 5 mins.

Step 4. Protein Removal

- a. Add 100 μ l B3 Buffer to the sample and vortex immediately for 10 seconds.
- b. Incubate on ice for 5 mins.
- c. Centrifuge at 14,000 x g for 3 minutes.
- d. Transfer the supernatant to a clear 1.5 ml microcentrifuge tube.

Step 5. DNA Binding

- a. Add 400 μ l BC Buffer to sample and shake vigorously.
- b. Place a BZ Column to a Collection Tube. Transfer the sample mixture to BZ Column.
- c. Centrifuge at full speed (13,000 rpm/17,900x g) for 30 seconds.
- d. Discard the flow-through and place BZ Column back in the Collection Tube.

Step 6. Wash

- a. Add 400 μ l W1 Buffer to BZ Column.
- b. Centrifuge at full speed (13,000 rpm/17,900x g) for 30 seconds. Discard the flow-through and place BZ Column back in the Collection Tube.
- c. Add 600 μ l W2 Buffer (ethanol added) to BZ Column.
- d. Centrifuge at full speed (13,000 rpm/17,900x g) for 30 seconds. Discard the flow-through and place BZ Column back in the Collection Tube.

Step 7. Dry

- a. Centrifuge at full speed (13,000 rpm/17,900x g) for 2 minutes to dry the column.

Step 8. Elution

- a. Place BZ Column to a clean 1.5 ml microcentrifuge tube (not provided).
- b. Add 50-90 μ l of pre-heated Elution Buffer (60 °C) into the center of the column matrix.
- c. Stand at 60 °C for 5 minutes.
- d. Centrifuge at full speed (13,000 rpm/17,900x g) for 2 minutes to elute purified DNA. Store the DNA fragment at 4 °C or -20 °C.

II. For Urine :

Step 1. Sample preparation

- a. Centrifuge the urine sample at 4000 x g for 5 minutes.
- b. Remove the supernatant completely and resuspend the pellet in 900 μ l B1 Buffer by pipetting.

Step 2. Cell Lysis

- a. Centrifuge 5 mins at 4000 x g. Remove the supernatant completely.
- b. Resuspend the cell with **50 µl B1 Buffer** by pipetting the pellet.
- c. **Add 300 µl B2 Buffer** to the sample and mix thoroughly by vortexing.
- d. Incubate at 60 °C for 10 mins until the sample lysate is clear. Invert the tube every 3 minutes during incubation.
- e. Preheat required Elution Buffer (50-90 µl per sample) in a 60 °C water bath.
- f. Centrifuge 5 mins at 3000 x g. Remove the supernatant completely.

Optional Step

If RNA-free genomic DNA is required, perform this optional step.

- a. Add 5 µl of RNase A (10 mg/ ml, not provided) to sample lysate and mix by vortexing.
- b. Incubate at room temperature for 5 mins.

Step 3. Protein Removal

- a. **Add 100 µl B3 Buffer** to the sample and vortex immediately for 10 seconds.
- b. Incubate on ice for 5 mins.
- c. Centrifuge at 14,000 x g for 3 minutes.
- d. Transfer the supernatant to a clear 1.5 ml microcentrifuge tube.

Step 4. DNA Binding

- a. **Add 400 µl BC Buffer to sample and shake vigorously.**
- b. Place a **BZ Column** to a Collection Tube. Transfer the sample mixture to BZ Column.
- c. Centrifuge at full speed (13,000 rpm/17,900x g) for 30 seconds.
- d. Discard the flow-through and place BZ Column back in the Collection Tube.

Step 5. Wash

- a. **Add 400 µl W1 Buffer** to BZ Column.
- b. Centrifuge at full speed (13,000 rpm/17,900x g) for 30 seconds. Discard the flow-through and place BZ Column back in the Collection Tube.
- c. **Add 600 µl W2 Buffer** (ethanol added) to BZ Column.
- d. Centrifuge at full speed (13,000 rpm/17,900x g) for 30 seconds. Discard the flow-through and place BZ Column back in the Collection Tube.

Step 6. Dry

- a. Centrifuge at full speed (13,000 rpm/17,900x g) for 2 minutes to dry the column.

Step 7. Elution

- a. Place BZ Column to a clean 1.5 ml microcentrifuge tube (not provided).
- b. **Add 50-90 µl of pre-heated Elution Buffer (60 °C)** into the center of the column matrix.
- c. Stand at **60 °C** for 5 minutes.
- d. Centrifuge at full speed (13,000 rpm/17,900x g) for 2 minutes to elute purified DNA.
Store the DNA fragment at 4 °C or -20 °C.

Product Use Limitation & Warranty

This product is intended to be used for life science research only.

It has not been approved for drug or diagnostic purpose.

YEASTERN' s products should not be resold, modified for resale, or used to manufacture commercial products without written approval by YEASTERN. YEASTERN guarantees the performance of all products in the manner described in our protocol. The purchaser must determine the suitability of the product for its particular use. Should any product fail to perform satisfactorily due to any reason other than misuse, YEASTERN will replace it free of charge.