



Genomic DNA Extraction Kit (Tissue)2.0

Product Use Limitation & Warranty

This product is intended to be used for life science research only. It has not been approved for drug or diagnostic purpose. YEASTERN's products should not be resold, modified for resale, or used to manufacture commercial products without written approval by YEASTERN. YEASTERN guarantees the performance of all products in the manner described in our protocol. The purchaser must determine the suitability of the product for its particular use. Should any product fail to perform satisfactorily due to any reason other than misuse, YEASTERN will replace it free of charge.

Ver. Q0125

No part of these protocols may be reproduced in any form or by any mean, transmitted, or translated into a machine language without the permission of YEASTERN BIOTECH CO., LTD.

Address: 6F-3, 23 Lane 169, Kang Ning St., Shijr, Taipei, 22180 Taiwan.
Tel: +886-2-2695-3922 **Fax:** +886-2-2695-3979
Email: yeastern@yeastern.com **Website:** www.yeastern.com

Copyright© 2017 All rights reserved. Yeastern Biotech Co., Ltd.
Copyright© 2017 All rights reserved. Yeastern Biotech Co., Ltd.

Cat. No.
SYG111-001
FYG111-100
FYG111-300

Genomic DNA Extraction Kit (Tissue)2.0

(SYG111-001/ FYG111-100/ FYG111-300) Store at RT/4 °C Ver. Q0125

Sample : 30 mg of fresh animal tissue
25 mg of paraffin-embedded tissue

Yield : Up to 50 µg

Contents

Item	SYG111-001 (4 preps)	FYG111-100 (100 preps)	FYG111-300 (300 preps)
T1 Buffer	1.5 ml	35 ml	95 ml
T2 Buffer	0.5 ml	12 ml	35 ml
W1 Buffer	2 ml	45 ml	125 ml
W2 Buffer	300 µl x 2	15 ml	25 ml x2
Elution Buffer	1 ml	10 ml	30 ml
GZ Column	4 pcs	100 pcs	300 pcs
Collection Tube	4 pcs	100 pcs	300 pcs
User Manual	1	1	1

Buffer Preparation

- Add ethanol (96-100%) to the W2 Buffer prior to first use.

Buffer	SYG111-001 (4 preps)	FYG111-100 (100 preps)	FYG111-300 (300 preps)
W2 Buffer	300 µl x 2	15 ml	25 ml x 2
Ethanol (96 - 100%)	1.2 ml x 2	60 ml	100 ml x 2

Important Notes

1. Buffers provided in this system contain irritants. Wear gloves and lab coat when handling these buffers.
2. Add ethanol (96- 100 %) to W2 Buffer when first open.
3. Prepare dry bath or water bath before the operation.
4. Resolve any precipitate by warming at 37°C.

Description

The Genomic DNA Extraction Kit (Tissue)2.0 is designed for rapid extraction of pure genomic DNA from fresh animal tissue or paraffin-embedded tissue. Efficiently remove cellular debris and inhibitors, this kit using column-type tube in purification process through three simple steps of binding, washing and then elution for the safe and convenient extraction of high-purity genomic DNA. The entire process can be completed in less than 1 hour without phenol/chloroform, and the final product can be used in PCR or other downstream experiments.

Additional Requirements

For General procedure

- a. Proteinase K (10 mg/ml)
- b. Absolute ethanol

For Paraffin-Embedded Tissue:

- a. Xylene

Purification Protocols

Step 1. Sample preparation

For Fresh Tissue

- a. Cut off 30 mg of fresh animal tissue.
- b. Grind the sample under liquid nitrogen to a fine powder with pestle and mortar.
- c. Transfer to a 1.5 ml tube.

For Paraffin-Embedded Tissue

- a. Slice up to 25 mg of paraffin-embedded tissue and transfer to a 1.5 ml tube.
- b. Add 1 ml xylene (not provided) and vortex vigorously and incubate at room temperature for 10 minutes.
- c. Vortex every 2 minutes during incubation.
- d. Centrifuge 3 minutes at 14,000 x g. Remove the supernatant.
- e. Add 1 ml absolute ethanol (not provided) and mix by vortexing.
- f. Centrifuge 3 minutes at 14,000 x g. Remove the supernatant.
- g. Repeat step (e-f) one more time.
- h. Open the tube and incubate at the 37 °C for 15 minutes.

Step 2. Cell Lysis

- a. Add 300 μ l of T1 Buffer and 20 μ l of proteinase K (10 mg/ml, not provided) to the sample.
- b. Incubate the sample at 60 °C for 30 minutes until the sample lysate is clear. Invert the tube every 5 minutes during incubation.
- c. Pre-heat the Elution Buffer at 75 °C.

Optional step:

- a. Add 5 μ l of RNase A (10 mg/ml, not provided) to sample lysate and mix by vortexing.
- b. Incubate at room temperature for 5 minutes.

If RNA-free genomic DNA is required, perform this optional step.

Step 3. Protein Removal

- a. Add 100 μ l T2 Buffer to the sample and shake vigorously.
- b. Centrifuge 14,000 x g for 3 minutes.
- c. Add 300 μ l absolute ethanol and shake vigorously.

Step 4. DNA Binding

- a. Place GZ Column with a Collection Tube.
- b. Transfer the sample mixture to the GZ Column and centrifuge at 14,000 x g for 30 seconds.
- c. Discard the flow-through and place the GZ Column back in the Collection Tube.

Step 5. Wash

- a. Add 400 μ l of W1 Buffer to the GZ Column.
- b. Centrifuge at 14,000 x g for 30 seconds.
- c. Discard the flow-through and place the GZ Column back in the Collection Tube.
- d. Add 600 μ l of W2 Buffer (ethanol added) to the GZ Column.
- e. Centrifuge at 14,000 x g for 30 seconds.
- f. Discard the flow-through and return the GZ Column back to the Collection Tube.

Note: Make sure that ethanol (96-100%) has been added into W2 Buffer when first use.

Step 6. Dry column

- a. Centrifuge at 14,000 x g for 3 minutes to dry the GZ column.

Step 7. Elution

- a. Combine the GZ Column with a 1.5 ml eppendorf (not provided).
- b. Add 50-90 μ l of preheated Elution Buffer (75 °C) into the center of the column matrix.
- c. Stand 75 °C for 3 minutes.
- d. Centrifuge at 14,000 x g for 2 minutes to elute purified DNA.

Step 7. Store DNA

- a. Store the DNA fragment at 4°C or -20°C.

Product Use Limitation & Warranty

This product is intended to be used for life science research only. It has not been approved for drug or diagnostic purpose. YEASTERN' s products should not be resold, modified for resale, or used to manufacture commercial products without written approval by YEASTERN. YEASTERN guarantees the performance of all products in the manner described in our protocol. The purchaser must determine the suitability of the product for its particular use. Should any product fail to perform satisfactorily due to any reason other than misuse, YEASTERN will replace it free of charge.