



## Genomic DNA Extraction kit (Cultured Cell) 2.0

### Product Use Limitation & Warranty

This product is intended to be used for life science research only. It has not been approved for drug or diagnostic purpose. YEASTERN's products should not be resold, modified for resale, or used to manufacture commercial products without written approval by YEASTERN. YEASTERN guarantees the performance of all products in the manner described in our protocol. The purchaser must determine the suitability of the product for its particular use. Should any product fail to perform satisfactorily due to any reason other than misuse, YEASTERN will replace it free of charge.

Ver. Q0307

No part of these protocols may be reproduced in any form or by any mean, transmitted, or translated into a machine language without the permission of YEASTERN BIOTECH CO., LTD.

**Address:** 6F-3, 23 Lane 169, Kang Ning St., Shijr, Taipei, 22180 Taiwan.  
**Tel:** +886-2-2695-3922 **Fax:** +886-2-2695-3979  
**Email:** [yeastern@yeastern.com](mailto:yeastern@yeastern.com) **Website:** [www.yeastern.com](http://www.yeastern.com)

Copyright© 2017 All rights reserved. Yeastern Biotech Co., Ltd.  
Copyright© 2017 All rights reserved. Yeastern Biotech Co., Ltd.

**Cat. No.**  
**SYG117-001**  
**FYG117-100**  
**FYG117-300**

# Genomic DNA Extraction kit (Cultured Cell) 2.0

SYG117-001/FYG117-100/FYG117-300

Store at RT/4 °C Ver. Q0307

Sample : 10<sup>7</sup> Cultured Cell

Yield: Up to 50 µg

## Contents

	SYG117-001	FYG117-100	FYG117-300
N1 Buffer	4 ml	100 ml	100 ml x3
N2 Buffer	1.5 ml	35 ml	95 ml
N3 Buffer	2 ml	45 ml	125 ml
W1 Buffer	2 ml	45 ml	125 ml
W2 Buffer	300 ul x2	15 ml	25 ml x2
Elution Buffer	1 ml	10 ml	30 ml
GZ Column	4 pcs	100 pcs	300 pcs
Collection Tube	4 pcs	100 pcs	300 pcs
User Manual	1	1	1

## Buffer Preparation

- Add ethanol (96-100%) to the W2 Buffer prior to first use

	SYG117-001	FYG117-100	FYG117-300
W2 Buffer	300 ul x2	15 ml	25 ml x2
Ethanol (96 ~ 100%)	1.2 ml x2	60 ml	100 ml x2

## Additional Requirements

1. Ethanol (96-100%)
2. RNase A(Optional)
3. Microcentrifuge tubes

## Important Notes

1. Buffers provided in this system contain irritants. Wear gloves and lab coat when handling these buffers.
2. Add ethanol (96- 100 %) to W2 Buffer when first open.
3. Prepare dry bath or water bath before the operation.
4. Resolve any precipitate by warming at 37°C.

## Purification Protocol :

### Step 1. Sample Preparation

- a. Transfer up to  $10^7$  cultured cells to a microcentrifuge tube (not provided).
- b. Centrifuge 1 minute at 6,000 x g. Remove the supernatant completely.
- c. Resuspend the cell with 50  $\mu$ l N1 Buffer by pipetting.

### Step 2. Cell Lysis

- a. Add 300  $\mu$ l N2 Buffer to the sample and mix thoroughly by vortexing.
- b. Incubate at 60 °C for 10 mins until the sample lysate is clear.
- c. Invert the tube every 3 minutes during incubation.

#### **Optional Step.**

*If RNA-free genomic DNA is required, perform this optional step.*

*Add 5  $\mu$ l of RNase A (10 mg/ ml, not provided) to sample lysate and mix by vortexing. Incubate at room temperature for 5 mins.*

### Step 3. Protein Removal

- a. Add 400  $\mu$ l N3 Buffer to the sample and shake vigorously.
- b. Centrifuge 1 minute at 12,000 x g.

#### Step 4. DNA Binding

- a. Place a GZ column in Collection Tube.
- b. Transfer the supernatant to GZ column and centrifuge 30 seconds at 14,000 x g.
- c. Discard the flow-through and place GZ Column back in the Collection Tube.

#### Step 5. Washing

- a. Add **400 µl of W1 Buffer** into the **GZ Column**.
- b. Centrifuge at full speed (14,000 x g) for 1 minute.
- c. Discard the flow-through and place the **GZ Column** back in the **Collection Tube**.
- d. Add **600 µl W2 Buffer** (Ethanol added) in the **GZ Column**.
- e. Centrifuge at full speed (14,000 x g) for 1 minute.
- f. Discard the flow-through and place the **GZ Column** back in the **Collection Tube**.

#### Step 6. Dry column

- a. Centrifuge at full speed (14,000 x g) for 3 minutes to dry the column.

#### Step 7. Elution

- a. Place **GZ Column** to a clean 1.5 ml microcentrifuge tube.
- b. Add **50-90 µl of preheated Elution Buffer (70°C)** into the center of the column matrix.
- c. Stand at room temperature for 3 minutes.
- d. Centrifuge at full speed (14,000 x g) for 2 minutes to elute purified DNA.

## Step 8. Store DNA

- a. Store the DNA fragment at 4°C or -20°C.

## Product Use Limitation & Warranty

This product is intended to be used for life science research only. It has not been approved for drug or diagnostic purpose. YEASTERN' s products should not be resold, modified for resale, or used to manufacture commercial products without written approval by YEASTERN. YEASTERN guarantees the performance of all products in the manner described in our protocol. The purchaser must determine the suitability of the product for its particular use. Should any product fail to perform satisfactorily due to any reason other than misuse, YEASTERN will replace it free of charge.