



## HiYield Gel/PCR DNA Fragments Extraction Kit 2.0

### Product Use Limitation & Warranty

This product is intended to be used for life science research only. It has not been approved for drug or diagnostic purpose. YEASTERN's products should not be resold, modified for resale, or used to manufacture commercial products without written approval by YEASTERN. YEASTERN guarantees the performance of all products in the manner described in our protocol. The purchaser must determine the suitability of the product for its particular use. Should any product fail to perform satisfactorily due to any reason other than misuse, YEASTERN will replace it free of charge.

Ver. Q0308

No part of these protocols may be reproduced in any form or by any mean, transmitted, or translated into a machine language without the permission of YEASTERN BIOTECH CO., LTD.

**Address:** 6F-3, 23 Lane 169, Kang Ning St., Shijr, Taipei, 22180 Taiwan.  
**Tel:** +886-2-2695-3922 **Fax:** +886-2-2695-3979  
**Email:** yeastern@yeastern.com **Website:** www.yeastern.com

Copyright© 2017 All rights reserved. Yeastern Biotech Co., Ltd.  
Copyright© 2017 All rights reserved. Yeastern Biotech Co., Ltd.

**Cat. No.**  
**SYG206-001**  
**FYG206-100**  
**FYG206-300**

# HiYield Gel/PCR DNA Fragments Extraction Kit 2.0

(SYG206-001/FYG206-100/FYG206-300)

Ver. Q0308

Sample : 100  $\mu$ l PCR Product, 300 mg of Agarose Gel

Store at RT/4°C

Yield : Up to 50  $\mu$ g

## Contents

|                 | SYG206-001     | FYG206-100 | FYG206-300 |
|-----------------|----------------|------------|------------|
| EZ Buffer       | 2 ml           | 60 ml      | 80 ml x2   |
| W1 Buffer       | 2 ml           | 45 ml      | 125 ml     |
| W2 Buffer       | 300 $\mu$ l x2 | 15 ml      | 25 ml x2   |
| Elution Buffer  | 1 ml           | 10 ml      | 30 ml      |
| EZ Column       | 4 pcs          | 100 pcs    | 300 pcs    |
| Collection Tube | 4 pcs          | 100 pcs    | 300 pcs    |
| User Manual     | 1              | 1          | 1          |

## Buffer Preparation

- Add ethanol (96-100%) to the W2 Buffer prior to first use

|                     | SYG206-001     | FYG206-100 | FYG206-300 |
|---------------------|----------------|------------|------------|
| W2 Buffer           | 300 $\mu$ l x2 | 15 ml      | 25 ml x2   |
| Ethanol (96 ~ 100%) | 1.2 ml x2      | 60 ml      | 100 ml x2  |

## Important Notes

1. Buffers provided in this system contain irritants. Wear gloves and lab coat when handling these buffers.
2. Add ethanol (96- 100 %) to W2 Buffer when first open.
3. Prepare dry bath or water bath before the operation.
4. Resolve any precipitate by warming at 37°C.

# Purification Protocols

## I. Gel Extraction :

### Step 1. Gel Dissociating

- a. Excise the agarose gel slice containing relevant DNA Fragments.
- b. Transfer up to 300 mg (**do not over 300 mg**) of the gel slice into a microcentrifuge tube (not provided).
- c. **Add 500  $\mu$ l EZ Buffer** to the sample and mix by vortexing.
- d. Incubate at 60 °C for 10 minutes until the gel slice has been completely dissolved. During incubation, invert the tube every 2-3 minutes.
- e. Cool down the dissolved sample mixture to room temperature slowly.

### Step 2. DNA Binding

- a. Place a **EZ Column** in a Collection Tube.
- b. Apply 800  $\mu$ l of the sample mixture into the EZ Column.
- c. Centrifuge at 14,000 x g for 30 seconds.
- d. Discard the flow-through and place the EZ Column back in the Collection Tube.

**Note: If the sample mixture is more than 800  $\mu$ l, repeat this DNA Binding Step.**

### Step 3-1. Wash

- a. **Add 400  $\mu$ l W1 Buffer** to EZ Column.
- b. Centrifuge at 14,000 x g for 30 seconds.
- c. Discard the flow-through and place the EZ Column back into the Collection Tube.

### Step 3-2. Wash

- a. **Add 600  $\mu$ l W2 Buffer** (ethanol added) into the EZ Column.
- b. Centrifuge at 14,000 x g for 30 seconds. Discard the flow-through and place EZ Column back in the Collection Tube.

### Step 4. Dry

- a. Centrifuge again for 2 minutes at 14,000 x g to dry the column matrix.

### Step 5. Elution

- a. Transfer dried EZ Column into a new microcentrifuge tube (not provided).
- b. **Add 50 - 90 µl Elution Buffer** into the center of the column matrix. ( If DNA is larger than 5 kb, use preheated 60 °C Elution Buffer to improve the elution efficiency. )
- c. Stand for 2 minutes until Elution Buffer is absorbed by the matrix.
- d. Centrifuge for 2 minutes at full speed to elute purified DNA.

### Step 6. Store DNA

- a. Store DNA at 4 °C or -20 °C.

## II. PCR Clean Up :

### Step 1. Sample preparation

- a. **Add 500 µl EZ Buffer** to 100 µl PCR product and mix by vortexing.

### Step 2. DNA Binding

- a. Place a **EZ Column** in a Collection Tube .
- b. Apply 800 µl of the sample mixture into the EZ Column.
- c. Centrifuge at 14,000 x g for 30 seconds.
- d. Discard the flow-through and place the EZ Column back in the Collection Tube.

**Note: If the sample mixture is more than 800 µl, repeat this DNA Binding Step.**

### Step 3-1. Wash

- a. **Add 400 µl W1 Buffer** to EZ Column.
- b. Centrifuge at 14,000 x g for 30 seconds.
- c. Discard the flow-through and place the EZ Column back into the Collection Tube.

### Step 3-2. Wash

- a. **Add 600 µl W2 Buffer** (ethanol added) into the EZ Column.
- b. Centrifuge at 14,000 x g for 30 seconds. Discard the flow-through and place EZ Column back in the Collection Tube.

#### Step 4. Dry

- a. Centrifuge again for 2 minutes at 14,000 x g to dry the column matrix.

#### Step 5. Elution

- a. Transfer dried EZ Column into a new microcentrifuge tube (not provided).
- b. **Add 50 - 90 µl Elution Buffer** into the center of the column matrix. (If DNA is larger than 5 kb, use preheated 60 °C Elution Buffer to improve the elution efficiency. )
- c. Stand for 2 minutes until Elution Buffer is absorbed by the matrix.
- d. Centrifuge for 2 minutes at full speed to elute purified DNA.

#### Step 6. Store DNA

- a. Store DNA at 4 °C or -20 °C.

### Product Use Limitation & Warranty

This product is intended to be used for life science research only.

It has not been approved for drug or diagnostic purpose.

YEASTERN' s products should not be resold, modified for resale, or used to manufacture commercial products without written approval by YEASTERN. YEASTERN guarantees the performance of all products in the manner described in our protocol. The purchaser must determine the suitability of the product for its particular use. Should any product fail to perform satisfactorily due to any reason other than misuse, YEASTERN will replace it free of charge.