



Total RNA Kit (Plant)2.0

Product Use Limitation & Warranty

This product is intended to be used for life science research only. It has not been approved for drug or diagnostic purpose. YEASTERN's products should not be resold, modified for resale, or used to manufacture commercial products without written approval by YEASTERN. YEASTERN guarantees the performance of all products in the manner described in our protocol. The purchaser must determine the suitability of the product for its particular use. Should any product fail to perform satisfactorily due to any reason other than misuse, YEASTERN will replace it free of charge.

Ver. Q0125

No part of these protocols may be reproduced in any form or by any mean, transmitted, or translated into a machine language without the permission of YEASTERN BIOTECH CO., LTD.

Address: 6F-3, 23 Lane 169, Kang Ning St., Shijr, Taipei, 22180 Taiwan.
Tel: +886-2-2695-3922 **Fax:** +886-2-2695-3979
Email: yeastern@yeastern.com **Website:** www.yeastern.com

Copyright© 2017 All rights reserved. Yeastern Biotech Co., Ltd.
Copyright© 2017 All rights reserved. Yeastern Biotech Co., Ltd.

Cat. No.
SYG308-001
FYG308-100
FYG308-300

Total RNA Kit (Plant)2.0

(SYG308-001/ FYG308-100/ FYG308-300) Store at RT/4 °C Ver. Q0125

Sample : 100 mg of Tissue

Yield : Up to 30 µg

25 mg of dry plant Tissue

Contents

Items	SYG308-001 (4 preps)	FYG308-100 (100 preps)	FYG308-300 (300 preps)
PR Buffer	4 ml	110 ml	105 ml x3
W1 Buffer	2 ml	45 ml	125 ml
W2 Buffer	300 ul x2	15 ml	25 ml x2
Elution Buffer	1 ml	10 ml	30 ml
RZ Column	4 pcs	100 pcs	300 pcs
Collection Tube	4 pcs	100 pcs	300 pcs
User Manual	1	1	1

Buffer Preparation

- Add ethanol (96-100%) to the W2 Buffer prior to first use.

Buffer	SYG308-001 (4 preps)	FYG308-100 (100 preps)	FYG308-300 (300 preps)
W2 Buffer	300 ul x2	15 ml	25 ml x2
ethanol (96 ~ 100%)	1.2 ml x2	60 ml	100 ml x2

Additional Requirements

1. β - Mercaptoethanol
2. RNase-free microcentrifuge tubes
3. Isopropanol
4. ethanol (96-100%)

Important Notes

1. Buffer contains chaotropic salt is harmful and irritant agent. Wear gloves and lab coat when handling these buffers.
2. Use sterile, RNase-free pipet tips and microcentrifuge tubes. Wear a lab coat and disposable gloves to prevent RNase contamination.
3. Make sure the starting sample amount is under the limit.
4. Add ethanol (96- 100 %) to W2 Buffer when prior to the initial use.
5. All purification steps should be carried out at room temperature.
6. All centrifugation should be carried out in a table-top microcentrifuge at $>12000 \times g$ (10,000-14,000 rpm, depending on the rotor type).

Purification Protocols

Step 1. Sample preparation

- a. Cut off 100 mg of fresh plant tissue or 25 mg of dry plant tissue.
- b. Grind the sample under liquid nitrogen to a fine powder by using a mortar and pestle.

Step 2. Cell Lysis

- a. Add 1 ml PR Buffer and 10 μ l of β - Mercaptoethanol to the sample in the mortar and grind the sample until it is completely dissolved.
- b. Transfer the sample mixture to a RNase-free microcentrifuge tube and incubate at 75°C for 30 minutes. (Invert the tube every 10 minutes.)
- c. Centrifuge at 2-8°C at 14,000 $\times g$ for 10 minutes and transfer the supernatant to a new microcentrifuge tube.
- d. Add 1/2 volume of isopropanol to the sample and shake vigorously.
- e. Place a RZ column in Collection Tube.

Step 3. RNA Binding

- a. Transfer the sample mixture (up to 700 μ l once) to RZ column and centrifuge 30 seconds at 14,000 $\times g$.
- b. Discard the flow-through and place RZ Column back in the Collection Tube.

Step 4. Wash

- a. Add 400 μ l W1 Buffer to RZ Column.
- b. Centrifuge at 14,000 x g for 30 seconds.
- c. Discard the flow-through and place RZ Column back in the Collection Tube.
- d. Add 600 μ l W2 Buffer (ethanol added) to RZ Column.
- e. Centrifuge at 14,000 x g for 30 seconds.
- f. Discard the flow-through and place RZ Column back in the Collection Tube.

Step 5. Dry column

- a. Centrifuge at 14,000 x g for 2 minutes to dry the RZ column.

Step 6. Elution

- a. Place RZ Column to a clean 1.5 ml microcentrifuge tube (not provided).
- b. Add 50-90 μ l of Elution Buffer into the center of the column matrix.
- c. Stand at room temperature for 2 minutes.
- d. Centrifuge at 14,000 x g for 2 minutes to elute purified RNA.

Step 7. Store RNA

- a. Store the RNA fragment at -80°C.

Product Use Limitation & Warranty

This product is intended to be used for life science research only. It has not been approved for drug or diagnostic purpose. YEASTERN' s products should not be resold, modified for resale, or used to manufacture commercial products without written approval by YEASTERN. YEASTERN guarantees the performance of all products in the manner described in our protocol. The purchaser must determine the suitability of the product for its particular use. Should any product fail to perform satisfactorily due to any reason other than misuse, YEASTERN will replace it free of charge.