Sharp Protein Markers III

Recommended storage conditions:

-20°C for one year Avoid multiple freeze/thaw cycles

Loading volume: Mini gel / blotting: 5 ul/well Large gel / blotting: 10 µl/well


Size range: 10~315 kDa

Fragments: 12 proteins Package size: 500 µl*1 vial

- The 72 kDa and 315 kDa reference bands are orange.
- The 52 kDa and 10 kDa reference bands are green.
- The protein marker is supplied in gel loading buffer and is ready-to-use. DO NOT BOIL, further dilution or addition of any reducing agent.
- After running the gel, stain the gel or perform Western blot transfer as desired



Sharp Protein Markers III

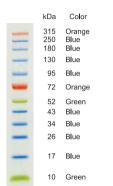

Mini gel loading volume: 5 µl/well 4~20% SDS-PAGE Gel



Yeastern Biotech Co., Ltd.