



Product Use Limitation & Warranty

This product is intended to be used for life science research only. It has not been approved for drug or diagnostic purpose. YEASTERN's products should not be resold, modified for resale, or used to manufacture commercial products without written approval by YEASTERN. YEASTERN guarantees the performance of all products in the manner described in our protocol. The purchaser must determine the suitability of the product for its particular use. Should any product fail to perform satisfactorily due to any reason other than misuse, YEASTERN will replace it free of charge.

Ver. R0426

No part of these protocols may be reproduced in any form or by any mean, transmitted, or translated into a machine language without the permission of YEASTERN BIOTECH CO., LTD.

Address: 6F-3, 23 Lane 169, Kang Ning St., Shijr, Taipei, 22180 Taiwan.
Tel: +886-2-2695-3922 **Fax:** +886-2-2695-3979
Email: yeastern@yeastern.com **Website:** www.yeastern.com

Copyright© 2017 All rights reserved. Yeastern Biotech Co., Ltd.
Copyright© 2017 All rights reserved. Yeastern Biotech Co., Ltd.

Total RNA Kit (Blood/ Bacteria)

Cat. No.
SYG306-001
FYG306-100
FYG306-300

Total RNA kit (Blood/ Bacteria)

(SYG306-001/FYG306-100/FYG306-300)

Store at RT/4 °C Ver. R0426

Sample : 10^9 x Bacteria Cells / 300 μ l of blood

Yield : Up to 30 μ g

Contents

	SYG306-001	FYG306-100	FYG306-300
BR1 Buffer	4 ml	110 ml	100 ml x 3
BR2 Buffer	2 ml	45 ml	125 ml
BR3 Buffer	1 ml	25 ml	65 ml
W1 Buffer	2 ml	45 ml	125 ml
W2 Buffer	300 μ l x2	15 ml	25 ml x2
Elution Buffer	1 ml	10 ml	30 ml
RZ Column	4 pcs	100 pcs	300 pcs
Collection Tube	4 pcs	100 pcs	300 pcs
User Manual	1	1	1

Add ethanol (96-100%) to the W2 Buffer prior to first use:

	SYG306-001	FYG306-100	FYG306-300
W2 Buffer	300 μ l x2	15 ml	25 ml x2
Ethanol (96 - 100%)	1.2 ml x2	60 ml	100 ml x2

Additional Requirements

1. β – Mercaptoethanol
2. Lysozyme Buffer (20 mg/ ml lysozyme)
3. RNase-free microcentrifuge tubes
4. 70% Ethanol

Important Notes

1. Buffer contains chaotropic salt is harmful and irritant agent.
2. Use sterile, RNase-free pipet tips and microcentrifuge tubes. Wear a lab coat and disposable gloves to prevent RNase contamination.
3. Make sure the starting sample amount is under the limit.
4. Add ethanol (96-100%) to W2 Buffer prior to the initial use.
5. All purification steps should be carried out at room temperature.
6. All centrifugations should be carried out in a table-top microcentrifuge at $>12000 \times g$ (10,000-14,000 rpm, depending on the rotor type).

Step 1. Sample preparation

I. For Gram- Bacteria :

- a. Transfer up to 10^9 bacteria cells to a microcentrifuge tube.
- b. Centrifuge at 12,000 xg for 1 min and remove the supernatant completely.
- c. **Add 200 μ l of BR3 Buffer** to the tube and resuspend the cell pellet by vortexing or pipetting.
- d. Incubate at room temperature for 5 minutes.
- e. **Add 300 μ l of BR2 Buffer** and 3 μ l β -Mercaptoethanol, mix by vortexing.
- f. Incubate at room temperature for 5 minutes.

II. For Gram+ Bacteria :

- a. Transfer up to 10^9 bacteria cells to a microcentrifuge tube.
- b. Centrifuge at 12,000 xg for 1 min and remove the supernatant completely.
- c. **Add 200 μ l of Lysozyme Buffer** (20 mg/ml lysozyme) to the tube and resuspend the cell pellet by vortexing or pipetting.
- d. Incubate at room temperature for 10 minutes.
- e. **Add 300 μ l of BR2 Buffer** and 3 μ l β -Mercaptoethanol, mix by vortexing.
- f. Incubate at room temperature for 5 minutes.

III. For Blood :

- a. Transfer up to 300 μ l of blood to a RNase-free microcentrifuge tube.
- b. **Add 900 μ l of BR1 Buffer**, then mix by inverting.
- c. Incubate the mixture on ice for 10 minutes, and invert every 5 minutes.
- d. Centrifuge at 4°C at 4,000 xg for 5 minutes.
- e. Remove the supernatant completely.
- f. **Add 100 μ l of BR1 Buffer** and resuspend the pellet by pipetting.
- g. **Add 400 μ l of BR2 Buffer** and 4 μ l β -Mercaptoethanol, mix by vortexing.
- h. Incubate at room temperature for 5 minutes.

Step 2. Cell Lysis

- a. Centrifuge at 14,000 x g for 10 minutes.
- b. Transfer the supernatant to a clean microcentrifuge tube.

Step 3. RNA Binding

- a. **Add 500 μ l of 70% ethanol** and shake vigorously.
- b. Place a **RZ Column** in **Collection Tube**.
- c. Transfer the sample mixture (up to 700 μ l once) to **RZ Column** and centrifuge 1 minute at 14,000 xg.
- d. Discard the flow-through and place **RZ Column** back in the **Collection Tube**.

Step 4-1. Washing

- a. **Add 400 μ l W1 Buffer** to **RZ Column**.
- b. Centrifuge at 14,000 x g for 30 seconds.
- c. Discard the flow-through and place **RZ Column** back in the **Collection Tube**.

Step 4-2. Washing

- a. Add 600 µl of W2 Buffer (ethanol added) into the column.
- b. Centrifuge at 14,000 x g for 30 seconds.
- c. Discard the flow-through and place the RZ Column back in the Collection Tube.

Step 5. Drying

- a. Centrifuge at 14,000 rpm for 2 minutes to dry the column matrix.

Step 6. RNA Elution

- a. Place RZ Column to a clean 1.5 ml microcentrifuge tube.
- b. Add 50-90 µl of preheated Elution Buffer (75°C) into the center of the column matrix.

Step 7. Purify RNA

- a. Stand at room temperature for 3 minutes.
- b. Centrifuge at 14,000 xg for 2 minutes to elute purified RNA.

Step 8. Store RNA

- a. Store the RNA fragment at -80°C.

Product Use Limitation & Warranty

This product is intended to be used for life science research only. It has not been approved for drug or diagnostic purpose. YEASTERN' s products should not be resold, modified for resale, or used to manufacture commercial products without written approval by YEASTERN. YEASTERN guarantees the performance of all products in the manner described in our protocol. The purchaser must determine the suitability of the product for its particular use. Should any product fail to perform satisfactorily due to any reason other than misuse, YEASTERN will replace it free of charge.